D

Anleitung CHECKIT[®] Nitrat

AF 525

15 52 50

0,20 - 20 mg/l Nitrat - Stickstoff (N)

0,88 - 88 mg/l Nitrat (NO₃)

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GB

Instructions for using the CHECKIT® Nitrate

AF 525

15 52 50

0.20 - 20 mg/l as N

0.88 - 88 mg/l as NO₃

1. Fill the Nitrate Test tube with the water sample under test to the mark (20 ml).
2. Add one level spoonful of Nitrate Test Powder and one Nitrate Test Tablet. Do not crush the tablet.
3. Replace the screw cap and shake well for one minute, then allow to stand for about 1 minute.
4. Gently invert the tube 3 or 4 times, afterwards allowing a 2 minute standing period for final settlement.
5. Carefully decant the clear water into the left-hand compartment of the CHECKIT® if results as N are required, or the right-hand compartment if results as NO₃ are required. Fill to the 10 ml mark.
6. The centre compartment should be filled with untreated sample to compensate for any inherent colour or turbidity present in the sample.
7. Add to the 10 ml of treated sample one Nitrite LR Tablet, crush with a clean stirring rod and replace the stopper.
8. Hold the stopper firmly in place and repeatedly invert the CHECKIT® until the tablet is fully dissolved. Allow to stand for 10 minutes.
9. After this standing period give the CHECKIT® a final shake and compare the colour produced against the standards, using daylight, selecting the nearest colour match.
10. The range of the test may be altered by simply changing the volume of the sample, and making up with distilled or deionised water as indicated below:

Important

1. Contents of nitrite in the sample react as well and can lead to false test results.
2. The test is not valid if the colour produced in the water sample is outside the range of the scale.
3. The center compartment is for compensation of any cloudiness and colour in the water sample.

Direct Reading			2 to 10 mg/l as N		Test Range:			
using a 20 ml sample			8.8 to 44 mg/l as NO ₃		4 to 20 mg/l as N			
			N mg/l NO ₃		17.6 to 88 mg/l as NO ₃			
N	mg/l	NO ₃	N	mg/l	NO ₃	N	mg/l	NO ₃
			10		44			
1.0		4.40	8		35.2	20		88
0.8		3.52				16		70.4
0.6		2.64	6		26.4	12		52.8
0.4		1.76	4		17.6	8		35.2
0.2		0.88	2		8.8	4		17.6

Test Range:

Simply follow steps 1 to 9 of the instructions above. Results are as displayed on the front of the Comparator, reproduced above.

Add 2 ml of sample to Nitrate-Test tube and make up to 20 ml volume with distilled or deionised water. Follow steps 2 to 9 of instructions above. Revised test results are as shown above.

Add 1 ml of sample to Nitrate-Test tube and make up to 20 ml volume with distilled or deionised water. Follow steps 2 to 9 of instructions above. Revised test results are as shown above.

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