## D

### Anleitung CHECKIT® Phosphat

AF 540

15 54 00

0,5 - 4,0 mg/l von PO<sub>4</sub>

0,16 - 1,30 mg/l von P

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## GB

### Instruction for using the CHECKIT<sup>®</sup> Phosphate (low range)

AF 540

15 54 00

0.5 - 4.0 mg/l as PO<sub>4</sub>

0.16 - 1.30 mg/ ass P

1. Remove the lid (see picture 1), rinse the 3 compartments of the CHECKIT<sup>®</sup> with the water sample under test and fill both outer compartments to the 10 ml mark. The centre compartment is also filled to compensate for any inherent colour or turbidity present in the sample.
2. For measuring the concentration of Phosphate as PO<sub>4</sub> use the right-hand compartment. For measuring the concentration of Phosphate as P use the left-hand compartment.
3. Add one Phosphate No.1 LR Tablett (see picture 2) to the relevant compartment, crush with a clean stirring rod and mix to dissolve. Then add a Phosphate No.2 Tablett to the same compartment and again crush and mix to dissolve.
4. Replace the lid (see picture 3) and allow to stand for 10 minutes for full colour development.
5. After this standing period, give the CHECKIT<sup>®</sup> a final shake and compare the colour produced against the standards using daylight, selecting the nearest colour match.
6. Read the Phosphate concentration in mg/l as PO<sub>4</sub> or P alongside the corresponding standard.
7. Clean the CHECKIT<sup>®</sup> thoroughly after each test.

### Important

1. The water sample under test should be below 25°C.
2. Boiling water should be filtered before testing.
3. Cloudiness and colour in the water sample can be influenced by the acidic reaction. In this case add one DPD - Acidifying tablet to the water sample in the center compartment. Proceed to test as above.
4. The test is not valid if the colour produced in the water sample is outside the range of the scale.
5. The center compartment is for compensation of any cloudiness and colour in the water sample.
6. Conversion factors:  
 $\text{mg/l as P}_2\text{O}_5 = \text{mg/l PO}_4 \times 0.75$   
 $\text{= mg/l P} \quad \times 2.29$

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