Thermostatschränke Thermostatically controlled incubators Etuves réfrigérées Armarios termostaticos Estufas reguladas por termóstato

AL654/654G ; AL655/655G
AL656/656G ; AL658/658G

DE	Betriebsanleitung	2 – 13
GB	Instruction Manual	14 – 25
FR	Mode d'emploi	26 – 37
ES	Instrucciones	38 – 49
PT	Manual de instruções	50 – 61

Inhalt

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Operating instructions

Content

1. Brief description	15
2. Function	15
3. Initial operation	15
3.1 Unpacking the appliance	15
3.2 Environmental protection	15
3.3 Mounting procedures	16
3.4 Installation of the appliance	17
3.5 Electrical connections	17
3.6 Drill holes	18
4. Adjustment and operation	19
4.1 Operating controls and indicators	19
4.2 Setting temperature values	20
4.3 Detecting a power failure	21
4.4 Temperature control and accuracy	21
4.5 Other operating controls	22
5. Maintenance and service	23
5.1 Cleaning	23
5.2 Shut down / non-use	23
5.3 Preparation for transport	23
6. Troubleshooting and repair	23
7. Technical data	24

1. Brief description

The thermostat-controlled incubators are designed to allow continuous regulation of temperature for various fields of application, e.g.

- 20 °C BOD₅ measurement
- 4 °C storage of sewage samples
- 25 °C enzyme activity (TTC-test)
- 37 °C colony count

All temperature control problems in the usual range of 2 °C to 40 °C are solved by the factory set universal plug-in controller which is maintenance-free and reliable. This appliance is not suitable for storing flammable liquids, acids and lyes.

2. Function

The inner temperature of the fully insulated incubator is controlled exactly by an intergrated corrosion-proof temperature probe which connects or switches-off either the cooling unit or the heating element.

The measured inner temperature is directly indicated by a display. The temperature range of 2 °C to 40 °C can be set in increments of 0,1 °C with the aid of two tactile response keys. The keyboard is protected by a rugged foil. Aircirculation is provided by 2 axial flow fans with a discharge width of 160 cm² and an output of 300 m³/h.

3. Initial operation

3.1. Unpacking the appliance

Inspect the appliance immediatly after unpacking for any damage which might have occurred during shipment. If any items are missing or damaged please contact the carrier and file a claim for damages. Contents of the delivery should be checked against the delivery note.

3.2. Environmental protection information

Please dispose of the packaging correctly! Take it to the nearest official disposal point instead of throwing it in the garbage can! Take proper care that the reffridgerant circuit pipes of the worn-out appliance are not damaged before it is taken to a suitable environmentally recognised refuse dump. This will ensure that the coolant contained within the appliance is not released into the environment but recycled.

3.3. Mounting

Conversion of the right-hand door to a left-hand door

- The appliance is usually delivered with the door mounted right-hand.
- Close the door and remove the hinge bracket ❶
- Remove the door in a downward movement
- Remove the upper hinge pin ❷ and insert it on the opposite side


Only AL654 / AL654G / AL655G / AL656G / AL658G:

- Hang the door on its hinge pin and close the door
- Mount the hinge bracket ❶ on the left side of the incubator (having first removed the plastic cover).
- Remove the handle ❸ from the left side of the door
- Remove the plastic cover from the right side of the door and screw the door handle on

Only AL655 / AL656 / AL658:

- Rotate the door 180°, transfer door handle ❹ and locking latch ❺ on the door
- Remove plastic cover ❻
- Hang the door on its hinge pin and close the door
- Mount the hinge bracket ❶ by screwing it to the left side of the incubator

Check the fit of the door, adjust it, if necessary, using the oblong slots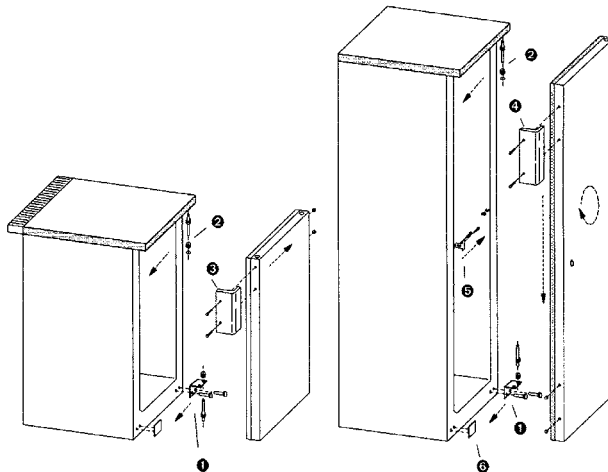

AL654 / AL654G / AL655G / AL656G / AL658G

AL655 / AL656 / AL658

3.4. Installation of the appliance

When choosing a location for the appliance the following points should be taken into account:

- no direct exposure to sunlight
- no proximity to any form of heating
- well ventilated rooms

The floor of the working area should be flat and level. Any unevenness can be compensated with the aid of the two adjusting feet (wrench size 22 mm).

The rear side of the appliance may be placed directly against the wall. The convection grids on the rear side of the appliance are not to be covered (e.g. table tops etc).

The carrying capacity of the retractable grids are about 45 kg.

Overall dimensions of the appliances in mm:

	Height	Width	Depth
AL654	850	600	600
AL654G	885	600	600
AL655	1216	600	600
AL655G	1216	600	600
AL656	1590	600	600
AL656G	1590	600	600
AL658	1516	752	710
AL658G	1516	752	710

3.5. Electrical connections

Connect the power cord to a shock-proof wall outlet. Any repositioning of the wall outlet may only be undertaken by a trained specialist. The VDE recommendations must be observed.

The socket must be protected with a fuse of at least 10 amps, which must not be in the rear section of the unit and must be easily accessible.

	Line voltage	Power consumption
AL654	230V/50Hz	200 VA
AL654G	230V/50Hz	210 VA
AL655	230V/50Hz	230 VA
AL655G	230V/50Hz	240 VA
AL656	230V/50Hz	230 VA
AL656G	230V/50Hz	240 VA
AL658	230V/50Hz	230 VA
AL658G	230V/50Hz	240 VA

Prior to putting the thermostat-controlled incubator into operation we recommend a thorough inner and outer cleaning of the appliance (see paragraph "cleaning" on page 23).


3.6 Drill holes

It is possible to provide drill holes for our thermostatic cabinets. On request, the drill holes may be made by the Production Department for new orders of a Thermostatic Cabinet. Alternatively, the customer may drill the holes on-site at a later date.

Please note: **In both cases, drilled holes will void the guarantee on the accuracy of the temperature's stability!** So long as the procedure is carried out correctly, statutory rights will not be affected.

The dimensions and possible drill areas are described in the graphic and respective tables below. In addition, the following should be taken in to consideration:

- Maximum hole size available by our production team: Ø 30 mm
- Wall thickness: 35 mm
- Initially, each hole is drilled from the inside to the outside with a diameter of approx. 3 mm. Then the hole can be drilled from the outside to the desired diameter / size.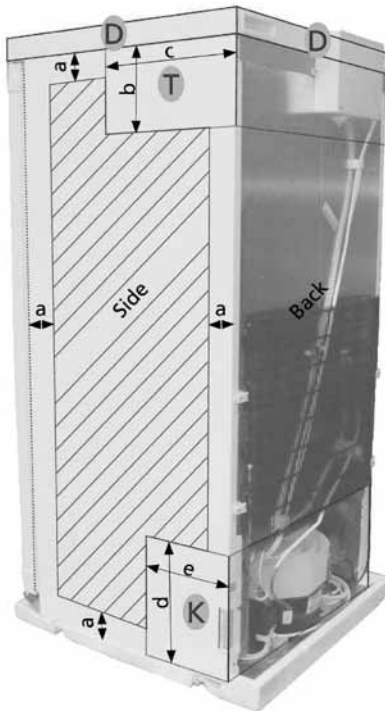

Available drilling area

a	External margins (mm)
All models	30

D	Cover height (mm)
AL654	70 (*)
Other models	80

(*) suitable for under bench applications, removable cover

T Area of Temperature Control Unit		
Model	b Height (mm)	c Depth (mm)
AL654(G)	230	330
AL655(G)	280	
AL656(G)	280	
AL658(G)	330	

K Area of Compressor Unit		
All models	d Height (mm)	e Depth (mm)
	230	230

The information provided is approximate. Slight variations may occur depending on the model.



4. Adjustment and operation

4.1. Operation controls and indicators

The unit is switched on and off via the On/Off switch.

The „Up” key is used to increase the temperature


The „Down” key is used to decrease the temperature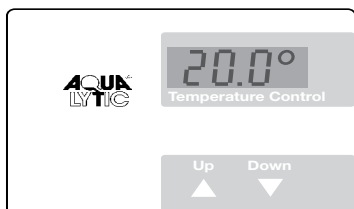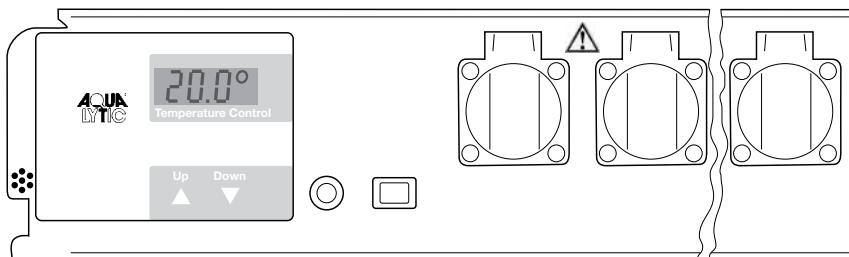

The illustration shows the actual temperature of 20.0 °C.

4.2. Setting the temperature values

When the unit is switched on, the display will initially show briefly the type of oven and a version number. It will then display the preset temperature, together with the symbol for the specified figure (3 horizontal lines). After 5 seconds, this will be replaced by the current, actual temperature.

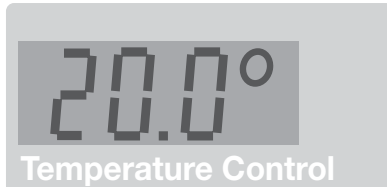The actual temperature can be read-off at any time (see also illustration of the display at the bottom of the previous page).

All temperatures between 2 °C and 40 °C can be set accordingly.

The unit is preset in the factory to a working temperature of 20° C. If this is required to be 25° C, for example:

Press the „Up“ key once. The display will show the specified figure which has been set, together with the symbol for the specified figure (3 horizontal lines). Now press the „Up“ key again, or hold it down until the display shows the new figure required (25° C in this example). Then release the key. The display will show the new setting for 5 seconds and will then display the actual, current temperature.

If a lower working temperature is required, the same procedure applies, using the „Down“ key. The temperature can be adjusted in steps of 0.1° C between 2° C and 40°C.

If you wish to read the specified temperature but without changing it, simply press the „Up“ or „Down“ key once and release it, so that the display shows the specified figure instead of the current temperature. The specified figure cannot be changed unless one of the keys is pressed while the display is already showing the specified figure.

The preset temperature will not be lost if there is a power failure or if the unit is switched off. The data are stored in an internal electronic memory and are called up automatically when power is restored or the unit is switched on.

Trouble shooting

If the verification of the temperature accuracy is not successful the following should be taken under consideration:

- The environmental temperature during the verification has to be between +10 and +32°C (climate class).
- A certain circulation of air is necessary for the compressor to work properly.
- The incubator should not be opened too often and / or for longer period of time.
- Check the function of the two fans by visual inspection.
- Check the cooling device by setting 10°C and wait for a significant change of the temperature in the sample towards 10°C.
- Check the heating device by setting 40°C and wait for a significant change of the temperature in the sample towards 40°C.
- Depending on the sample temperature it can take up to 7 hours for temperature stability at the set-value.

Important:

The air temperature in the cabinet may fluctuate by several degrees centigrade. Nevertheless the actual temperature of the sample fluctuates only by a maximum of $\pm 0.5^{\circ}\text{C}$.

4.5 Other operation controls

On the front side of the temperature control unit there are covered shock-proof sockets for the connection of BOD's appliances or stirring apparatuses. The total load on the sockets must not exceed 5 amps.

The thermostatically controlled oven and laboratory sockets are protected by a fuse (6.3 A T) which is accessible from the front panel of the unit.

The control unit is connected by means of a 4-pole plug on the bottom of the incubator.

5. Maintenance and controls

5.1 Cleaning

Every 6 - 8 months the inside of the unit should be freed of dust. At the same time, retractable grids at the rear side of the appliance should be freed of dust using a dry dusting brush. Care must be taken that no cables are pulled out or tubing bent or damaged in the process.

For a thorough cleaning of the inside of the appliance (e.g. after a long period of storage) use lukewarm water containing some detergent or other mild domestic cleaner on the plastic surfaces.

5.2 Shut down/non-use

If the thermostat-controlled incubator remains unused over a long period it should be switched off (disconnect plug from the mains or switch off or remove the fuse) and cleaned (see above!). To avoid bad odours the door should remain open during this period of non-use.

Prior to all work on the inside of the unit the mains plug of the thermostat-controlled incubator must be pulled out of the mains connection (not just a switching-off at the mains).

The appliance has an automatic defrosting device. Defrosted water collects in an evaporator tray and is evaporated automatically. Attention must be paid that the defrosted water can drain off freely through the drainage opening in the back panel of the incubator. The drainage opening should, when necessary be cleaned with the aid of a suitable device (e.g. bottle brush).

5.3 Preparation for transport

The appliance must be sent shock-proof-packed (preferably in its original packing). The packing is to be labelled as correspondingly ("keep dry/attention, fragile").

6. Troubleshooting and repair

Attention !!!

Before undertaking any repair work disconnect plug from mains! Only trained specialists should work on electrical devices.

Although switched ON - no display or control function

Cause: safety fuse defect

Remove defect fuse and replace it with a new fuse (slow-blow 6,3 A).

Correct display of inside temperature but no regulation to the temperature setpoint

Cause: compressor of the cooling system or control unit is defect

Contact your dealer or supplier, if necessary send in the control unit or thermostat-controlled incubator for repair (original packing if possible!).

7. Technical data

Construction:	fully insulated incubator with universal plug-in temperature
Control unit operation:	robust, water protected foil front panel, 2 keys with tactile acknowledgement
Range of adjustment:	2 °C to 40 °C; adjustable in 0.1 °C-increments
Constancy of temperature:	20 °C, ± 1.0° C (see 4.4)
Climate class:	10 °C to 32° C
Ambient temperature:	10° - 35°C
Display:	3 digit for actual and target temperature, resolution 0.1 °C
Refrigerating unit:	compressor-system
Heating unit:	heating element
Fan:	axial, discharge capacity 300 m ³ /h
Power requirement:	230V/50Hz
Coolant:	R134a

AL654 (4 sockets)

Net capacity:	135 Ltr.
Power consumption:	200 VA 1,35 kWh / 24h (ambient temperature 25 °C, temperature temperature 20 °C)
Overall dimensions:	850 H x 600 W x 600 D mm with worktop 820 H x 600 W x 600 D mm without worktop
Inside dimensions:	700 H x 515 W x 430 D mm (with 3 retractable grids and 1 bottom grid)
Weight:	39.0 kg
Order Code:	42 82 00

AL654G (4 sockets)

Net capacity:	140 Ltr.
Overall dimensions:	885 H x 600 W x 600 D mm
Inside dimensions:	700 H x 515 W x 443 D mm
Power consumption:	210 VA ; 1,77 kWh / 24h (ambient temperatur 25 °C, target temperature 20 °C with interior lighting switched on (15 W))
Weight:	50.0 kg
Order Code:	42 82 10

same as AL654, door, however, double glazed insulated in anodized aluminium frame

AL655 (5 sockets)

Net capacity:	195 Ltr.
Power consumption:	230 VA 1,39 kWh / 24h (ambient temperature 25 °C, temperature temperature 20 °C)
Overall dimensions:	1216 H x 600 W x 600 D mm
Inside dimensions:	1045 H x 515 W x 415 D mm (with 4 retractable grids and 1 bottom grid)
Weight:	49,0 kg
Order Code:	42 82 20

AL655G (5 sockets)

Net capacity: 195 Ltr.
Power consumption: 240 VA ; 1,74 kWh / 24h (ambient temperature 25 °C, target temperature 20 °C with interior lighting switched on (15 W))

Weight: 66.5 kg
Order Code: 42 82 25

same as AL655, door, however, double glazed insulated in anodized aluminium frame

AL656 (6 sockets)

Net capacity: 280 Ltr.
Power consumption: 230 VA
1,20 kWh / 24h (ambient temperature 25 °C, temperature 20 °C)

Overall dimensions: 1590 H x 600 W x 600 D mm
Inside dimensions: 1450 H x 515 W x 415 D mm
(with 5 retractable grids and 1 bottom grid)

Weight: 64,5 kg
Order Code: 42 82 30

AL656G (6 sockets)

Net capacity: 280 Ltr.
Power consumption: 240 VA ; 2,05 kWh / 24h (ambient temperature 25 °C, target temperature 20 °C with interior lighting switched on (15 W))

Weight: 82.0 kg
Order Code: 42 82 35

same as AL656, door, however, double glazed insulated in anodized aluminium frame

AL658 (8 sockets)

Net capacity: 395 Ltr.
Power consumption: 230 VA
1,59 kWh / 24h (ambient temperature 25 °C, temperature 20 °C)

Overall dimensions: 1516 H x 752 W x 710 D mm
Inside dimensions: 1300 H x 652 W x 515 D mm
(with 5 retractable grids and 1 bottom grid)

Weight: 79,5 kg
Order Code: 42 82 40

AL658G (8 sockets)

Net capacity: 395 Ltr.
Power consumption: 240 VA ; 1,97 kWh / 24h (ambient temperature 25 °C, target temperature 20 °C with interior lighting switched on (15 W))

Weight: 98.5 kg
Order Code: 42 82 45

same as AL658, door, however, double glazed insulated in anodized aluminium frame

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