LOVIBOND

MINIKIT

GB: Total Hardness, 1,4 - 56 °e
D: Gesamthärte, 1,0 - 48 °dH
F: Durete totale, 2,0 - 80 °f
I: Durezza totale
E: Durezza total
P: Durezza Total
NL: Totale hardheid
DK: Total hårdhed

20 - 800 mg/l CaCO₃

AF 445
41 44 50

GB

Introduction

This Minikit is for determination of Total Hardness (as CaCO₃) in water. The test range is 20-800 mg/l CaCO₃.

Test measurement Total Hardness, i.e. the total content of calcium and magnesium hardness.

Contents

dispenser bottle
calibrated sample container, 100 ml
calibrated sample container, 30 ml
T - Hardness tablets, 50 pieces in foil
operating manual in 8 languages

Instructions

1. Fill the dispenser bottle with the water sample to be tested. The dispenser bottle is used to allow a controlled addition of sample into the graduated tube.
2. Rinse the calibrated sample container with the water sample, leaving a few drops in the bottom.
3. Add one T - Harness - Test tablet and allow to disintegrate completely. A **BLUE** colour will be produced.
4. Using the dispenser bottle, gradually fill the tube with the sample until the colour changes from

BLUE to VIOLET.

Shake the tube gently during addition to ensure mixing and be careful not to overshoot the colour change.

5. The calcium hardness is given by reading graduation corresponding to the water level in the tube.

Cleaning

Thoroughly rinse out the dispenser bottle and the calibrated sample container after each use.

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