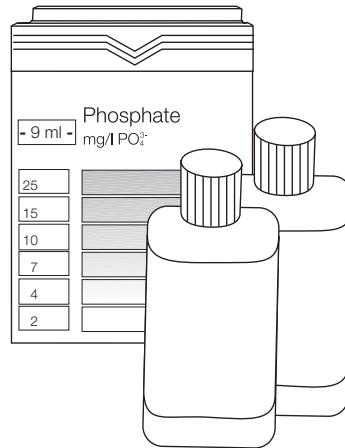
# Ortho-Phosphate $\text{PO}_4^{3-}$

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Bestimmung in  
Kesselwasser  
Determination  
in boiler water

Technical changes without notice  
Printed in Germany 02/20

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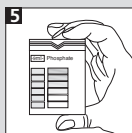
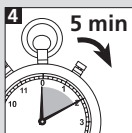
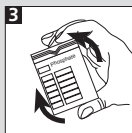
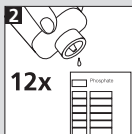
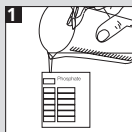
### Ortho-Phosphate $\text{PO}_4^{3-}$ (Phosphate-1/Phosphate-2)

- 1 Rinse comparator several times with the filtered water sample and fill the left and middle chamber of the comparator to the 9 ml mark.
- 2 Add 12 drops of reagent Phosphate-1 to the left chamber.
- 3 Replace the cap tightly and swirl gently. 2 Remove the cap, add 12 drops of reagent Phosphate-2. 3 Replace the cap tightly and swirl gently.
- 4 Wait for a colour reaction time of 5 minutes.
- 5 Match the colour fields against north daylight and read off the result as mg/l  $\text{PO}_4^{3-}$ .

### Notes

1. If the colour intensity of the sample is darker than the colour scale, the test has to be repeated with a diluted sample (for example 10 ml of sample + 10 ml of distilled water = dilution factor 2; the reading must be multiplied by 2).  
**Attention:** Dilute before adding reagent.
2.  $\text{SiO}_2$  concentrations higher 25 mg/l give increasing  $\text{PO}_4^{3-}$ -readings. Repeat test with a diluted sample (see point 1 above).
3. Citrates, oxalates, tartrates and similar agents reduce the sensitivity of the test.
4. Pyro-, meta-, and poly-phosphate do not react.  
 $1 \text{ mg/l } \text{PO}_4^{3-} \triangleq 0,75 \text{ mg/l } \text{P}_2\text{O}_5$ .