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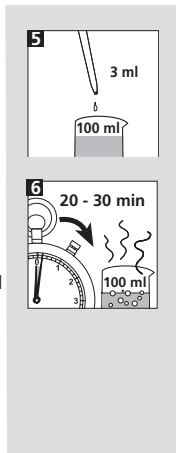
total-Phosphate

- 5 Add 3 ml concentrated hydrochloric acid to a 100 ml sample. Digest at 100°C for 20 to 30 minutes.
 6 After cooling down replace the evaporate water with deionized water. Continue with 1. The result is total-Phosphate in mg/l PO_4^{3-} . Reading B.

poly-Phosphate


Reading B - Reading A = poly-Phosphate in mg/l PO_4^{3-}

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


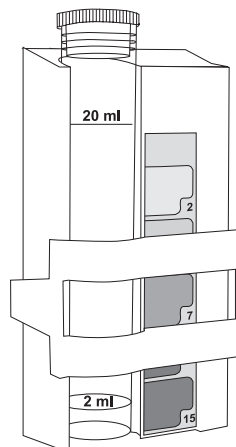
ortho-Phosphat (Phosphat-Reagenz I / Phosphat Reagenz II)

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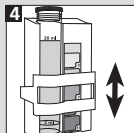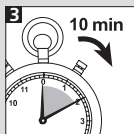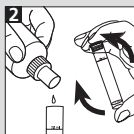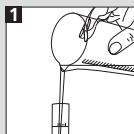Instructions for the determination of Phosphate

Instruments

Comparator unit Phosphate	Order Code 41 95 03
Test tube 2-20 ml	41 95 02

Reagents

Phosphate-Reagent I	Order Code (per 250 ml) 41 94 05
Phosphate-Reagent II	41 94 06

ortho-Phosphate (Phosphate-Reagenz I / Phosphate Reagenz II)

- 1 Fill a test tube to the 20 ml-mark.
- 2 Hold the dosage bottle vertical, remove the lid. By squeezing the bottle the top gets filled with 1 ml. The excess reagent returns into the bottle by loosening the pressure. Close the bottle immediately after use. Add 1 ml Phosphate-Reagent I in the test tube, close it with a lid. Swirl it gently. Add 1 ml Phosphate-Reagent II to the test tube, close it with a lid. Swirl it gently.
- 3 Wait 10 Minutes for full colour reaction.
- 4 Place the test tube in the comparator. Read off the result as mg/l PO_4^{3-} . Reading A.

Note

If the range of the comparator scale is exceeded, fill the test tube with the sample to the 10 ml / 5 ml / 2 ml-mark and fill up with deionized water to the 20 ml-mark. Continue with 2. Multiply the result with 2 at 10 ml / 4 at 5 ml / 10 at 2 ml.

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*) Bayer® AG, Leverkusen