CSB/COD-Reaktor

ET 125

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(GB) **Instruction Manual**
Page 8 - 12

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1. CSB-Reaktor ET 125

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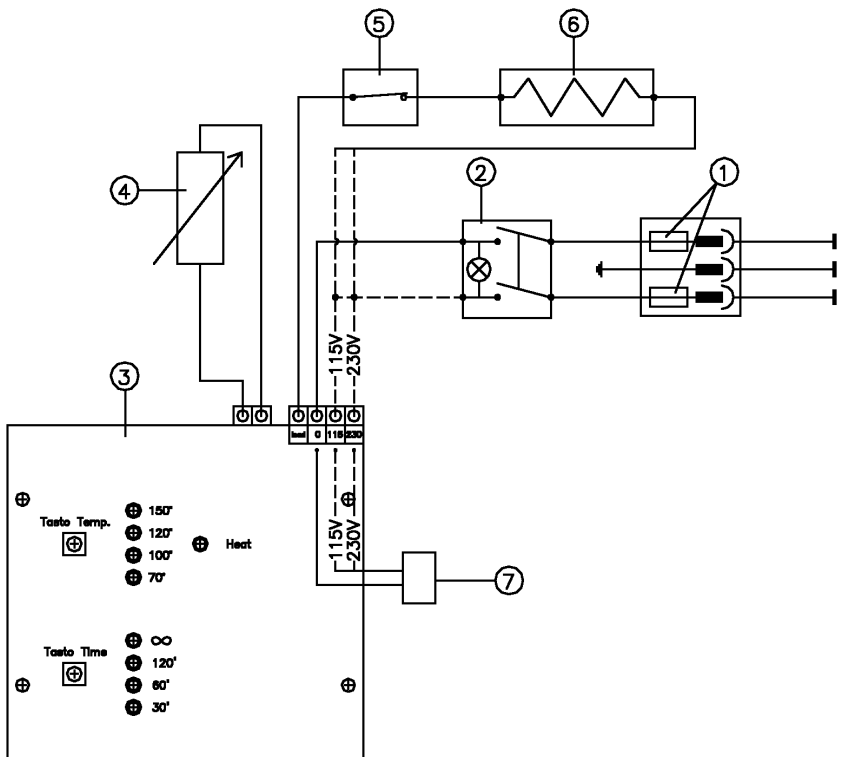
Contents

1. COD-Reactor ET 125

1.1. Important Notes	8
1.2. Safety Rules	8
1.3. Cleaning	8
1.4. Personal Protection Equipment	8
1.5. Measurement	9
1.6. Electric scheme	11

1. COD-Reactor ET 125

1.1. Important Notes

Read the information contained in the present manual before using the unit:

The manufacturer does not acknowledge any responsibility for an improper use of the equipment, not responding to directions for use.

The labels applied to the unit warn the user on the dangers to which he is exposed during the use or the maintenance. The labels must be left on the unit and substituted if they are no more readable.

Please use in all cases the protection cover, when heating
COD Reaction Cuvettes in Reactor.

Read all instructions before using the instrument.

1.2. Safety rules

1. The heating plate if programmed, may reach a temperature of 150 °C, this happens during the heating phase when the green LED is lighted, but also when it is not lighted during the cooling phase.
2. The materials used during the work must be compatible with the temperatures reached by the unit.

1.3. Cleaning

Always unplug the unit before cleaning. The heating plate must be cool. Use a damp cloth and not flammable, not aggressive detergents.

1.4. Personal Protection Equipment

The equipment used for personal protection must be compatible with the reached temperature and the dangers due to the working materials.

1. COD-Reactor ET 125

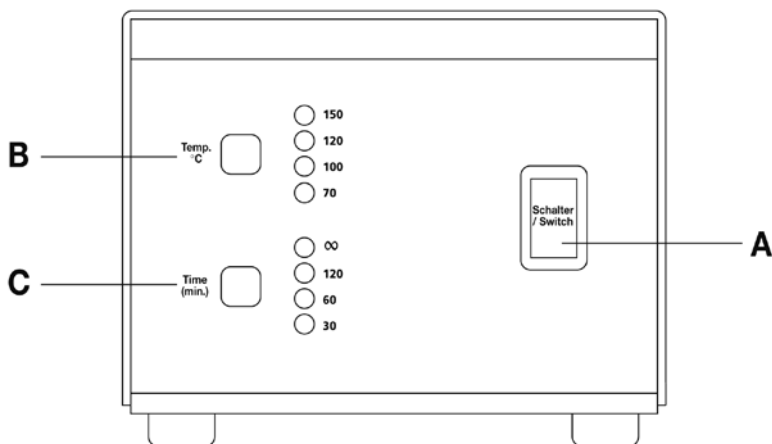
1.5. Measurement

1. Connect the instrument to power 230V/50 Hz (115V/60 Hz) by the cable and two pin plus ground.
2. The instrument is turned on by the mains switch (A) located on front panel. Automatically the instrument starts a work cycle with a temperature of 150 °C and a heating time of 120 minutes. The used values are shown by the lighting of corresponding LEDs on front panel. The heating time starts only when the reactor reaches the chosen temperature as shown by the lighting of the green LED (D) located on the upper part of front panel. The end of the heating period is warned by a beeper and the heating is turned off at the same time. For starting a new heating cycle it is necessary to turn off and on again the mains switch (A).
3. If a change in heating temperature or time is wanted, after turning on the instrument by the mains switch (A), the buttons (B) or (C) are to be pressed changing the values which are shown by the corresponding LEDs .

The possible values are:

Temperature 70/100/120/150 °C

Time 30 minutes or 60 minutes or 120 minutes or infinite
(no switching off = continuous operation at the chosen temperature)

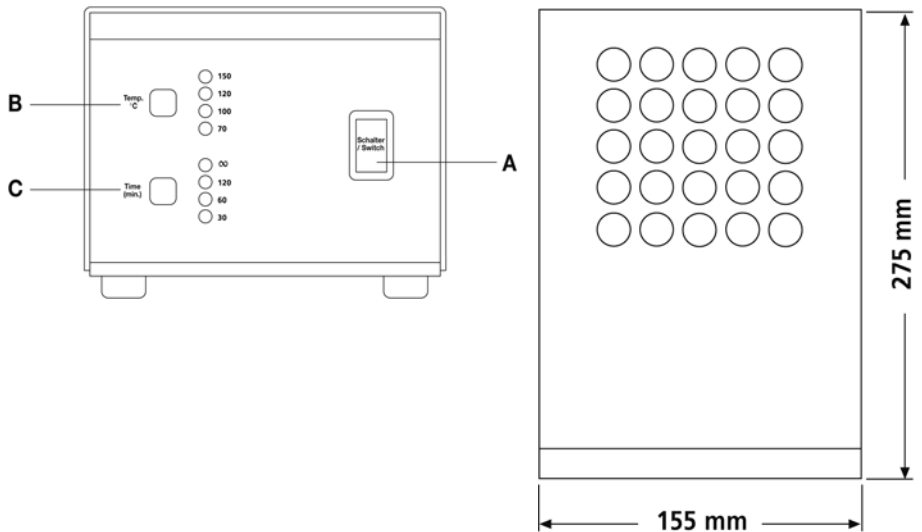
1. COD-Reactor ET 125

It is possible to change the values of temperature or time in any phase of the heating cycle. In this case the instrument starts again from zero and does not consider anymore the first choice.

4. The reactor is equipped by two retarded fuses (one spare) enclosed into the safety plug located on rear panel.
5. To clean the instrument we suggest to use tissue moistened with ethanol after disconnecting the plug from power.
6. The apparatus is planned and assembled according to safety standard IEC 1010 (1990) and modification 1 (1991).

Power	400 W
Weight	3,60 kg
Supply	230 V / 50 Hz
Supply	115 V / 60 Hz

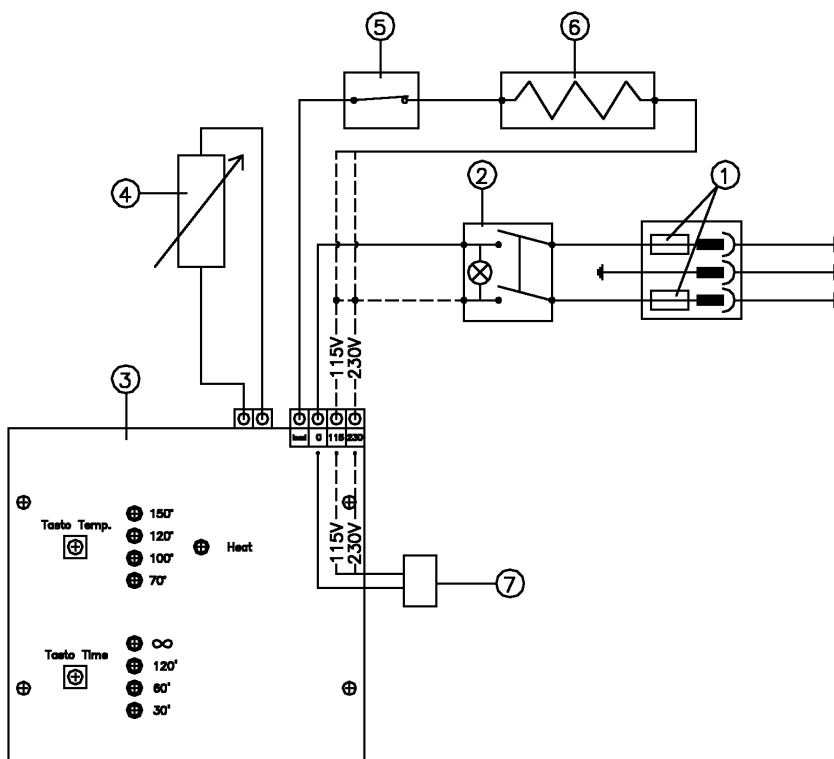
1. COD-Reactor ET 125

1.6. Electric scheme

Electric scheme 230 V/50 Hz (115 V/60 Hz)

1. 5 x 20 mm retarded fuse 3.15 A, (8 A)
2. Bipolar lighted switch
3. PT 100 probe
4. Resistance
5. Control Electronic Card
6. Security Thermostate

Technische Änderungen vorbehalten
Printed in Germany 06/09
We reserve the right to alter
or amend any of the items
contained herein without prior notice.

AQUALYTIC®
Postfach 41 02 53
44272 Dortmund
Germany
Telefon: (+49) (0)2 31/9 45 10 - 755
Telefax: (+49) (0)2 31/9 45 10 - 750